# LIEBHERR LOADING PLAN

### SCIENTIFIC REFRIGERATOR LKv 3910 / LKv 3913




## LOADING WITH PFIZER/BIONTECH TRAYS FOR 195 VIALS

#### **General Information**

The temperature-controlled Pfizer Cool Box enables the logistic of the Pfizer / BioNTech vaccine BNT162b2 at a temperature of -70°C by usage of dry-ice. One box/container can be filled with 5 trays, each tray with 195 vials. According to available information 5 vaccination doses can be taken from one vial.

The reusable Pfizer Cool Box consists of an outer carton and a payload sleeve for the five trays. The gap between outer carton and payload sleeve is filled up with dry-ice. With one dry-ice-loading the container keeps a temperature of -70°C for at least 10 days. The vaccines can be stored at +2°C to +8°C for max. 5 days before used.

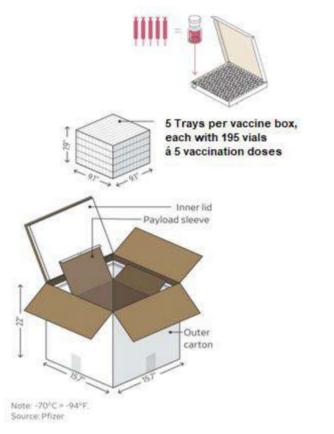


Source: https://si.wsj.net/public/resour ces/images/B3-HG530\_COLDVA\_300PX\_202009 03213118.jpg

Vaccine boxes / Payload (5 trays): height 7.9" = 20.1 cm x width 9.1" = 23.2 cm x depth 9.1" = 23.2 cmTransport container / outer carton: height 22" = 55.9 cm x width 15.7" = 39.9 cm x depth 15.7" = 39.9 cm



### LKv 3910 / LKv 3913 Loading Capacity Pfizer / BioNTech BNT162b2

**7 vaccine-boxes**, each vaccine box with 5 trays + **2 additional trays** (green marked in sketch), **overall 37 trays**; Each tray with 195 vials for 5 doses = **7.215 vials** / **36.075 doses** (or people that can get a vaccination) <u>Note:</u> for appliances with interior light 2 trays / 390 vials / 1.950 doses less (LKv 3913)

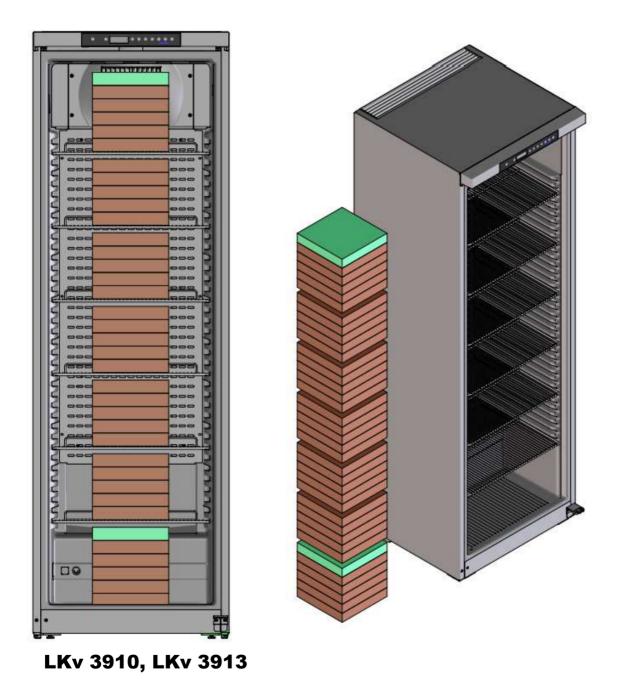

Disclaimer: All information regarding the Pfizer Cool Box as well as its dimensions are without guarantee. This information refers to existing press reports and may change due to further developments.



2020-11-24/HAU-PMB, lwlwar0